

### ISOLATION OF FECAL DNA

---



Vortex Bashing Beads Lysis  
Tube with fecal or soil sample

10,000 x g, 1 min



Filter lysate with Spin II  
A Filter

8000 x g, 1 min



Bind, Wash, Elute DNA  
with Spin Column

10,000 x g, 1 min



Filter DNA with Spin II  
B Filter

8000 x g, 3 min



Transfer eluted DNA

16,000 x g, 3 min



Isolated DNA

Technical Support

The troubleshooting guide is provided in the online protocol.  
For technical support please email us at [mbi.tech@meridianlifescience.com](mailto:mbi.tech@meridianlifescience.com) or visit [www.bioline.com](http://www.bioline.com)